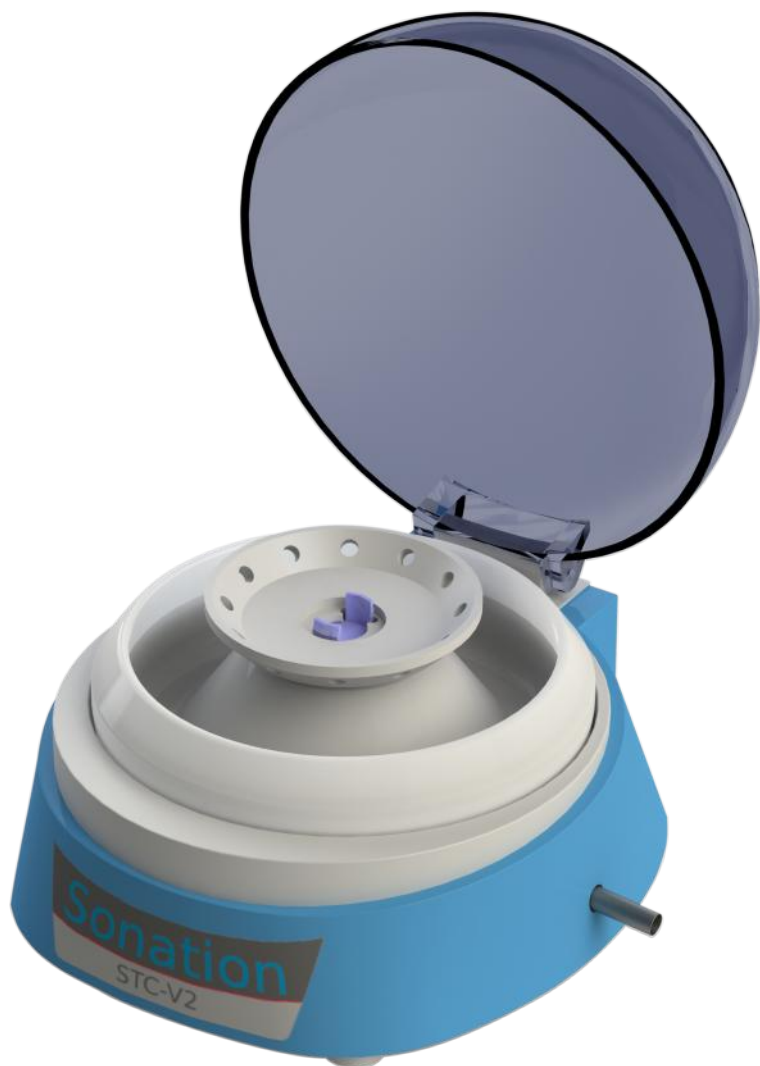
STC-V2 Zentrifuge für Stage Tipping STC-V2 Stage Tipping Centrifuge

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DESCRIPTION

Stage tipping is a sample preparation method developed at the Max Planck Institute for Biochemistry, department of proteomics and signal transduction in Munich - Germany. A video tutorial on the stage tipping method can be found under the following link:

<http://www.biochem.mpg.de/226863/Tutorials>

The Sonation Stage Tip Centrifuge is a modification of the TOMY Multi Spin lab centrifuge. The modification makes it possible to use open ended stage tips. The escaping solvent is collected and can be passed through a drain hose directly into a waste bottle.

This manual describes the handling and features of the modified centrifuge. For the basic operation of the centrifuge, please refer to the also included original users manual by TOMY.

OPERATION

Exchanging the Stage Tip Rotor

The Stage Tip Rotor can be unlocked and lifted by pressing the violet receptacle. In order to be able to centrifuge closed vials, also the original TOMY rotor can be used.

Inserting and removing the collecting funnel

The collecting funnel is held in place by magnets and the down-facing drain nipple. It can be easily taken out by lifting it up. When reinserting it must be ensured that it rests all around on the centrifuge base because only then it is ensured that no solvent can penetrate the housing. The drain nipple of the funnel locks into place with a noticeable "click" into the seal in the outlet of the centrifuge.

Mounting and changing the drain hose

The drain hose is simply pushed onto the drain tube. When laying the hose to the waste bottle care must be taken to ensure that the solvent can flow freely and that no back pressure forms. The hose must not be higher than the centrifuge outlet at any point. Otherwise, solvent may leak into the centrifuge or the collecting funnel may overflow. Because the hose is made of silicone, it can become brittle over time, depending on the solvents used. In this case, the hose should be replaced. Replacement hoses can be purchased from Sonation.

Inserting the stage tips and starting the centrifuge

The stage tips are inserted into the receiving holes of the rotor. If less than 12 stage tips are to be used, care must be taken that they are put in symmetrically in order to prevent imbalance during centrifuging.

If the centrifuge is connected to a wall outlet and the rear switch on, it is started by closing the lid and stopped again by opening the lid or switching the centrifuge off.

SPARE PARTS

STC-V2-CF	Collecting funnel with drain nozzle
STC-V2-STR	Stage Tip Rotor
STC-V2-DH	Drain hose